# Degassing





Fluidic Connections | Valves | Pumps | Degassers | Column Hardware | Manifolds

## **Degassing Solutions**



IDEX Health & Science degassers improve fluidic instrument precision and reliability by removing dissolved gases from fluids before they outgas and form problem-causing bubbles

We offer a broad portfolio of vacuum degassing assemblies to control bubbles in a wide range of system-fluids and flow-rates.

#### Degassing / Debubbling Membrane Materials

Widest Fluid Capability

Most Aqueous Based Capability



AF

AF based degassers offer the widest range of chemical compatibility and are best used with non-water-based system fluids.



Silicone

Silicone based degassers offer high flow rate capabilities for water-based systems to improve dispense accuracy and reliability.



#### Poridex

Poridex based debubblers provide rapid bubble removal for locations where bubble introduction cannot be avoided.

COMMON APPLICATIONS IMPROVED WITH DEGASSING

- Diagnostic Instrumentation: Clinical chemistry, immunoassay, hematology and molecular diagnostics
- Analytical Instrumentation: High performance liquid chromatography, sample handling and preparation
- Biotechnology: Next generation sequencing, sample handling, preparation and other microfluidic applications

### Importance of Degassing

Degassers avoid inaccurate sampling, maintain throughput and reduce errors by removing dissolved gases – even across complex fluid systems – before they outgas into bubbles. Utilizing degassing in your fluidic path is essential for consistent and accurate results.

Degassers greatly improve precision by degassing fluids far below the saturation point so instrument operations remain bubble-free. Without degassing, repeat sampling is required to detect inaccurate readings, which reduces throughput along your fluidic path. If bubble formation is not initially identified, sample assessment may be flawed. As the graph below demonstrates, degassers greatly increase accuracy for dispense volumes over time, in standard operation.



#### Dispense Volume Over 1,000 Dispenses\*

\* In most water-based applications

Temperature, pressure, reagent concentration and sock or agitation are all common variables that reduce the saturation point of systems fluids, resulting in outgassing that forms bubbles. IDEX Health & Science degassers manage efficiency across a wide range of flow rates and conditions with closed-loop vacuum control, to greatly enhance run-torun reliability and throughput, while maintaining dispense accuracy to avoid sample loss.



#### Solubility of Gas System Fluids

Any time system conditions shift in a way that reduces the gas solubility level of a fluid, outgassing occurs and bubbles form. If conditions change towards the lower right at any point in the fluid path, bubbles will be generated.

#### Learn More on Our Website

Explore our full list of degassing products, 360 degree interactive 3D images, complete specifications & more!

www.idex-hs.com/degassers



## **Degassing Methods**



### Degassers

Degassers eliminate bubbles before they form, which enhances instrument accuracy. Degassers offer the greatest advantage in complex systems, considering a single degasser can degas bulk fluid throughout an entire operation, regardless of the number of terminal flow paths.



The fluid stream enters the degassing chamber containing dissolved gas molecules and possibly bubbles

A vacuum is used to pull the dissolved gas molecules from the fluid stream

The fluid exiting the degasser is now free of dissolved gasses



#### OUTGASSING

Outgassing occurs when a solution's gas saturation point is altered due to the following causes:

- Capacity changes such as temperature increases or reduction in pressure
- Mixing different types of reagents or a change in reagent chemistry

#### Degassing in Bulk Fluids

Placing our degassers directly after the point which bulk fluids are being aspirated into an instrument will keep bubbles from forming throughout your system. Even small changes can result in trace amounts of outgassing to accumulate over time, in junctions and fittings. Bubbles introduced by loading and unloading bulk fluids, can be captured and removed by debubblers before causing system problems.





### Debubblers

Debubblers rapidly clear air pockets where bubble introduction cannot be avoided. Due to their small size, debubblers can be used along the fluidic path, after potential sources of bubble introduction, where they continuously remove bubbles, reduce downtime, and maintain optimal throughput.

As the fluid stream enters the debubbler, bubbles are trapped at the top of the chamber while fluid continues moving downstream

The vacuum source extracts the air bubbles and removes them from the fluid stream. An optional degassing tail can remove dissolved gasses in the fluid stream



The patented vertical design of IDEX Health & Science Debubblers trap bubbles at the top, removing them from the fluidic path.

#### Alternative to Flow Path Adjustment

Degassing replaces the need for manual flow path adjustments, and eliminates bubble build-up on probe tips – which cause level sensing errors – resulting in less reagent usage and increased throughput.

#### Debubblers for Optimal Throughput

Debubblers protect sensitive points of the fluid path to maintain throughput, which is the key objective of nearly all large diagnostic instruments. Bubbles are captured early, and purged to prevent erroneous repeated measurements.



#### BUBBLE INTRODUCTION

Bubble introduction occurs when air is brought into the flow path by the following causes:

- > Bulk fluid handling and changing
- > Sample injection
- > System leaks
- > Air permeable system tubing

## **Targeting Your Solution**

Complete Degassers DEGASSING CHAMBERS & VACUUM SYSTEMS





Degassers for

SILICONE MEMBRANE

Aqueous Based Fluids



Type of Degassing System Number of Independent Fluid Channels Dimensions (L x W x H)			Stand Alone Degassers	OEM Mini	OEM Mini Lite	250w Degassing Chamber	300w Degassing Chamber
			1 – 5	1 – 5	1-6	1	1
			10.4 x 2.9 x 5.2" (26.4 x 3.4 x 13.2 cm)	7.9 x 6.0 x 2.4" (20.0 x 15.2 x 6.1 cm)	6.8 x 2.1 x 4.0" (17.3 x 5.3 x 10.2 cm)	9.9 x 3.5 x 3.5" (25.1 x 8.9 x 8.9 cm)	3.8 x 3.5 x 3.5" (9.7 x 8.9 x 8.9 cm)
1	Where are the bubbles coming from?	Outgassing		~			✓
		Introduced Leaks		×			×
2	What type of fluid does your system use?	>50% Water Solution		•			~
		<50% Water Solution		<b>~</b>			×
3	What maximum flow rate is needed?			≤ 40 mL/min		≤ 250 mL/min	≤ 300 mL/min
4	Which vacuum system should be used?	Analytical Vacuum Control	Pump Included			×	
	Analytical for <10 mL/min Prep for >10 mL/min	Prep Vacuum Control				*	



For a full list of degassing products, please visit www.idex-hs.com/degassers

We are a strong force of committed people and innovative products for your fluidic pathway, continually increasing our product offering, expanding our market relevance by connecting to new customers, and positioning ourselves as global leaders in fluidics engineering.

## Worldwide Fluidics at IDEX Health & Science

As a global company, IDEX Health & Science has an international network of direct sales professionals and distribution partners in place to provide personal service to every customer. Our experts are ready to visit your operation, assess your fluidic needs, and develop solutions for your application challenges.

#### Corporate Responsibility

IDEX Health & Science is committed to preserving the environment. Our continuous improvement programs hold our facilities accountable to reduce waste, prevent pollution, and conserve resources. Many products comply with REACH and RoHS regulations.

#### North America

**IDEX Health & Science LLC** Bristol, CT, USA Carlsbad, CA, USA Middleboro, MA, USA Oak Harbor, WA, USA Rochester, NY, USA Rohnert Park, CA, USA Wallingford, CT, USA Asia ERC KK Saitama, Jap



#### For ordering and technical support, please contact:

 North America CustomerService.hs@idexcorp.com | +1 800 426 0191 | +1 360 679 2528

 Europe CustomerService.hsEurope@idexcorp.com | +49 1801 808 800 | +49 9377 9203-0

 Asia CustomerService.hsAsia@idexcorp.com | North Asia +86 10 6566 9090 | +86 21 5241 5599

 South Asia +91 22 66 43 55 50 | Southeast Asia & Oceania +65 6763 6633 | Japan +81 48 240 5750

 © 2015 IDEX Health & Science LLC. IDEX Health & Science LLC is a Unit of IDEX Corporation. | IDX2273

For More Information Visit idex-hs.com



Fluidic Connections | Valves | Pumps | Degassers | Column Hardware | Manifolds